

From Eye to Insight



Leica AirLab Icon Guide


Leica AirLab

- Download the Leica Airlab App from the Apple App store for iOS devices or from Google Play for Android devices. See note on next page before downloading!!




Leica AirLab

- If during the App download you get the below message regarding Leica AirLab accessing your Photos, please select OK. This simply allows the automatic creation of a Leica Collection in your photo Library and you will be able to share your AirLab images directly into the Leica Collection. If you select DON'T ALLOW, a Leica Collection will not be added and the mobile device operating system will not allow you to add a Leica Collection later.

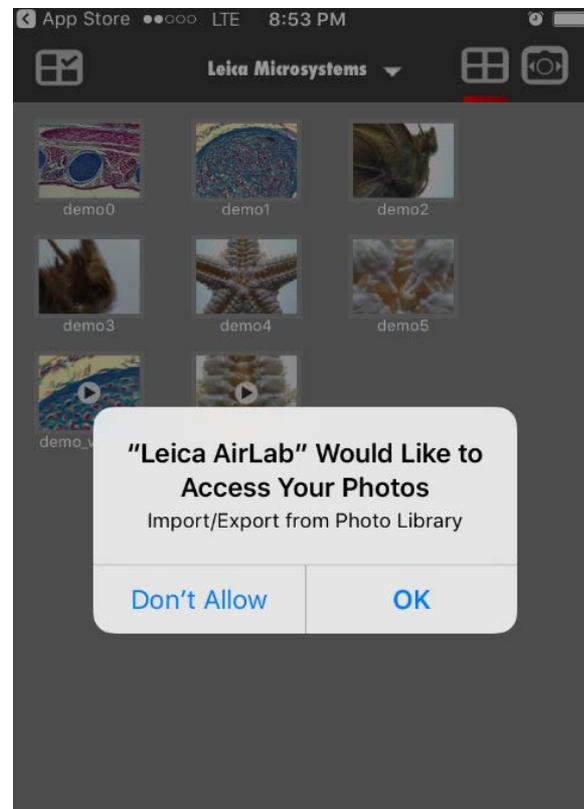


Leica AirLab

- Select your camera WiFi network in your device settings

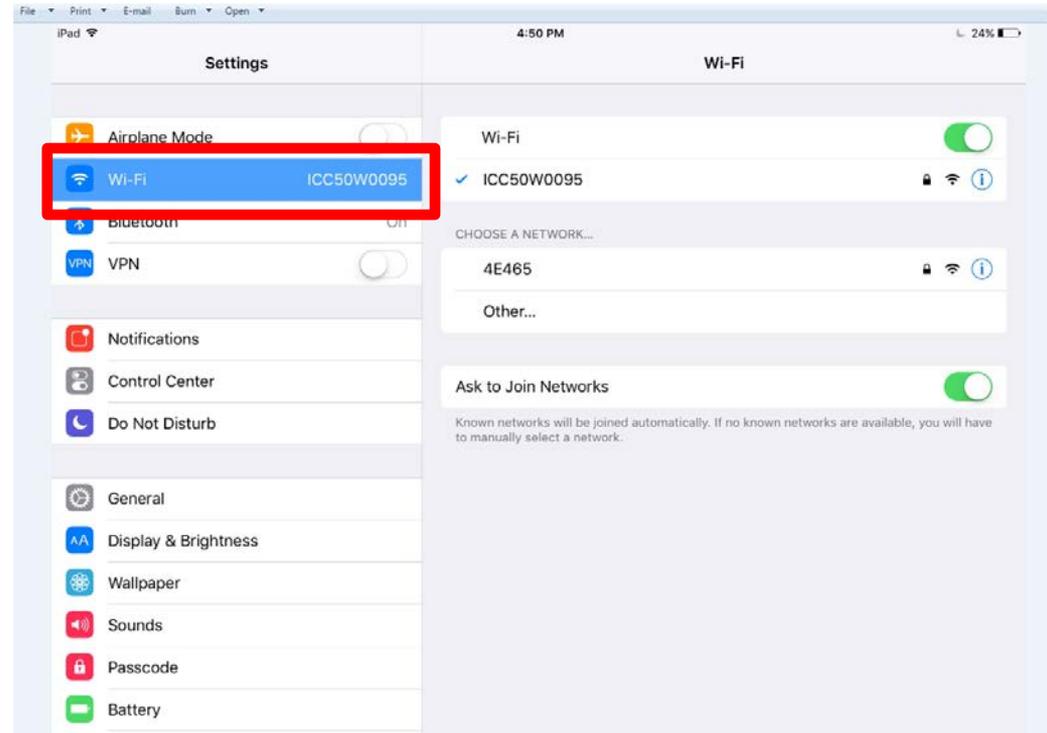

Leica AirLab

- Find AirLab in your App List of your device Settings and toggle on or off „Lite“ mode for allowing camera control or not

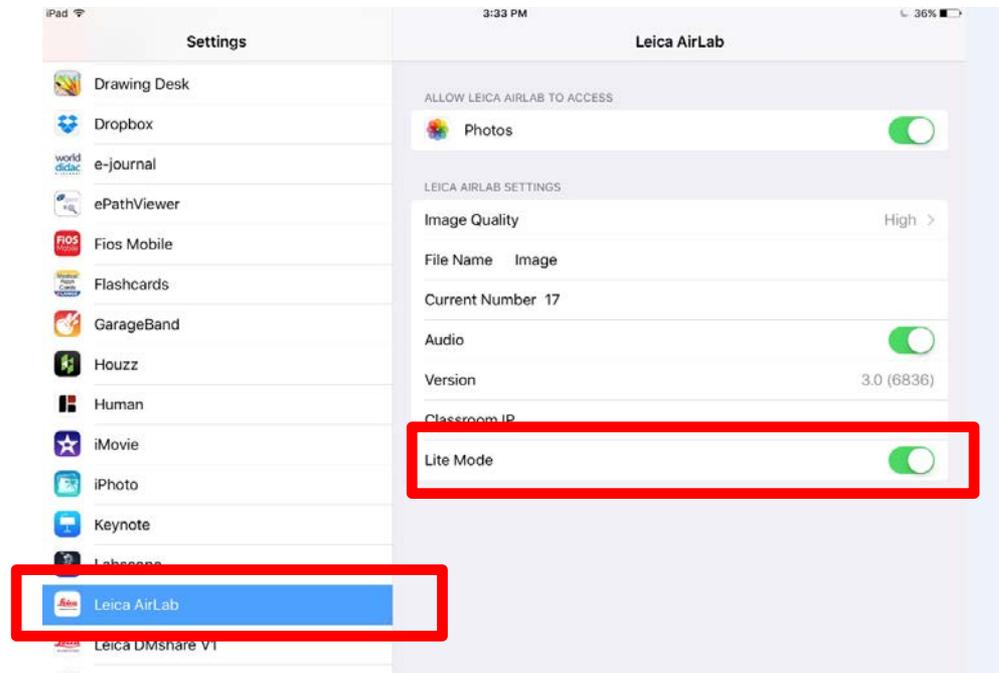



Leica AirLab

- Open the Leica AirLab App





Leica AirLab

- AirLab opens in Gallery View

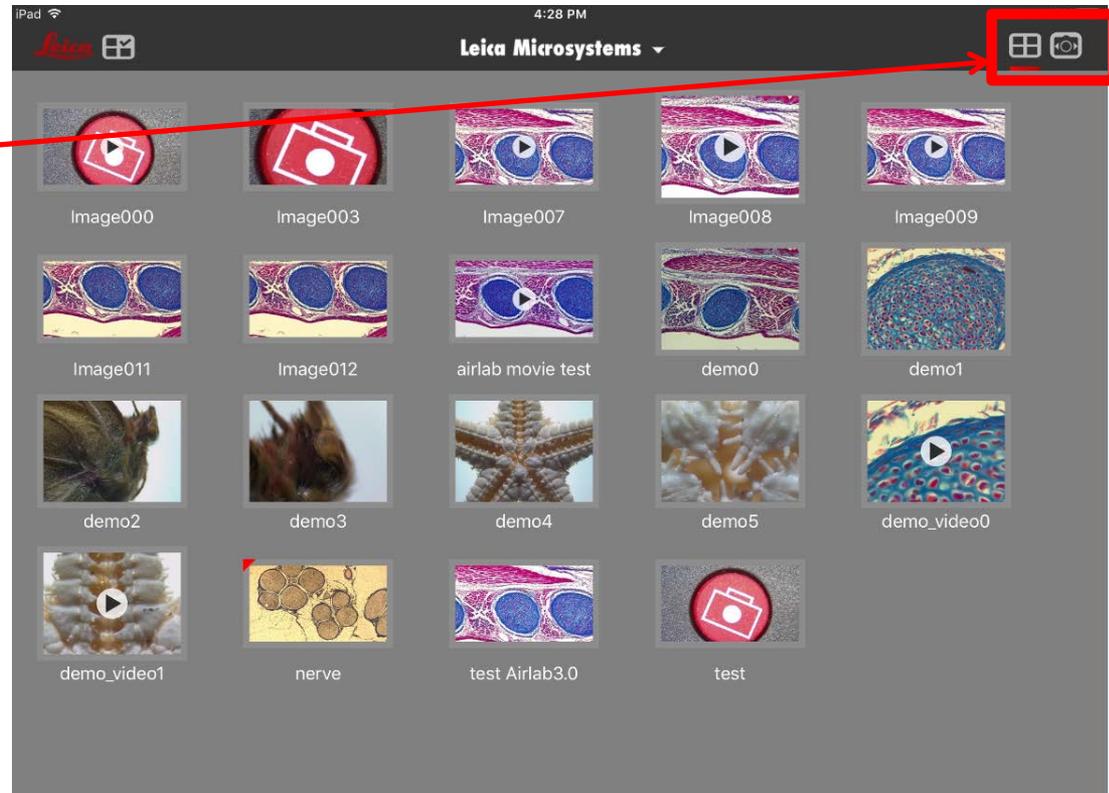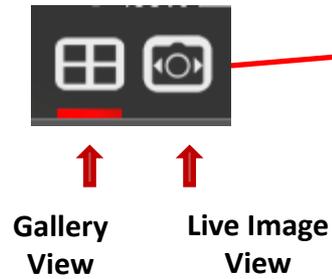

Collections

- Select/Create a Collection

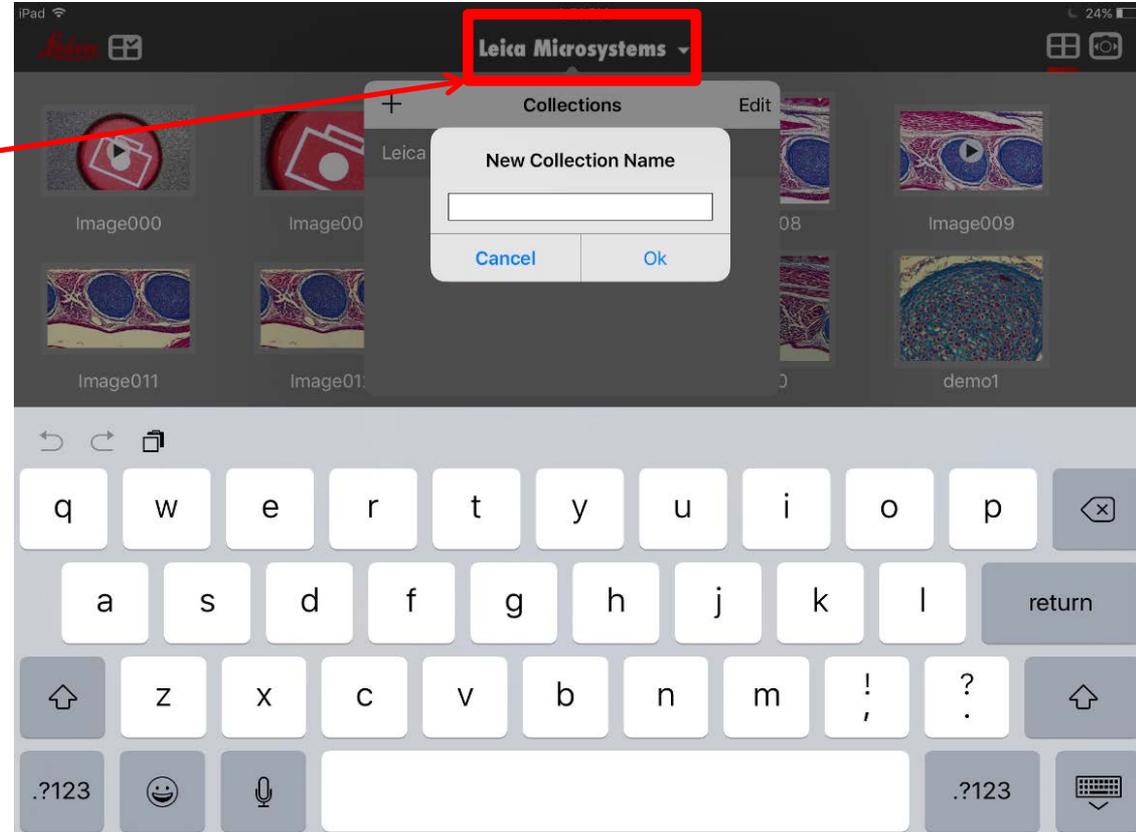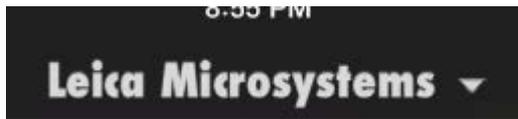

Select Live Image View



Add a camera by IP address/URL



Select Camera list view



Live Image View


- Select Live Image View

Tap an image to select it for full screen view.

Note that the thumbnails first appear as a Leica logo, then after a few seconds, the microscope image appears.



The thumbnails are updated approximately every 5 seconds.

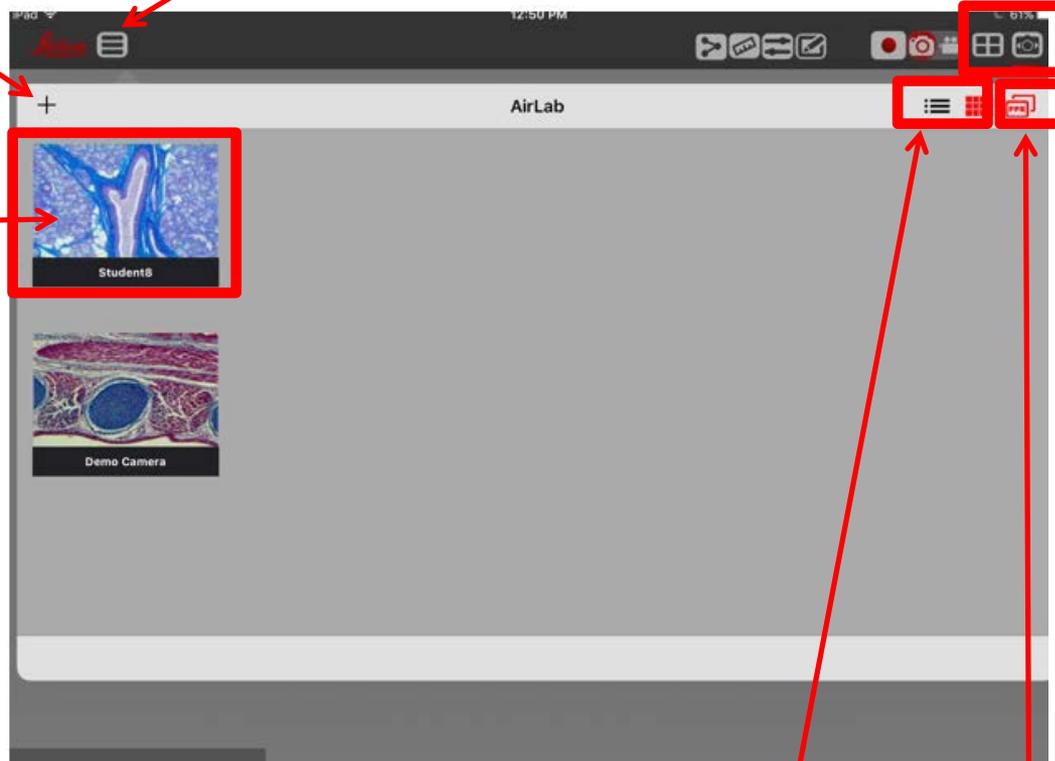

If you are connecting to a network and subnet where several cameras are connected, you will see thumbnails of all cameras connected to the network



Select thumbnail view or list view



Toggle on/off view of frame rate and live resolution

Calibrate



Calibration

- To calibrate the system select the calibration icon from the live image view
- A pop up shows default magnifications
- Tap the magnification you are using
- Tap „Set“
- The Live image is now Calibrated for the magnification you selected
- To Add a magnification, tap „Add: and another line will be added to the list
- To modify the Magnification, tap „Modify“



Calibrate



Calibration

- When you tap „Modify“ a line appears on the screen along with a pop up on the lower right
- Place an object of known dimension (ie: micrometer) under the microscope and stretch the line to the known dimension
- Tap on the pop up details to change the name, dimension, units of measure, and color
- Tap „OK“



Camera Settings

- Select the camera settings icon for control of the acquisition parameters



↑
Camera Settings



Camera Settings

- Camera Setting elements



Rotate /Flip the image

Auto exposure/Brightness
Manual exposure/Exposure time, Gain

Gamma, saturation, sharpening, auto white balance, region white balance, Live resolution

Return to original settings



Note when changing Live resolutions, the capture resolution will change to offer choices in the same aspect ratio

Capture Formats	Live Formats
Live	0.5MP (4:3)
720p (16:9)	0.8MP (4:3)
2.0MP (4:3)	720p (16:9)
2.5MP (4:3)	2.0MP (4:3)
1080p (16:9)	1080p (16:9)
5.0MP (4:3)	

Annotations and Measurements

- Select the annotation icon to get access to the toolbox



Annotations

Select all annotations



Mark and delete items




Drop downs for line, measurement, arrow, angle



Drop downs for circle, radius, ellipse





Drop downs for rectangle, free hand, text, insert image from your device photo library

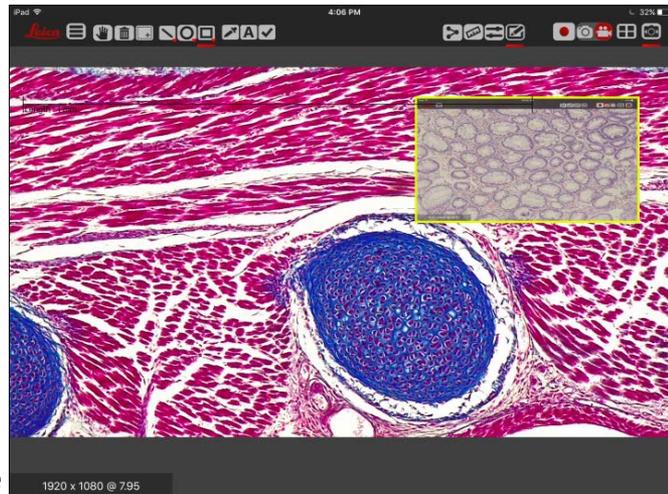

Annotations and Measurements

- Adjust font, color and size of annotations
- Turn on or off annotations, scale bar, measurement values, # of decimals for measurements

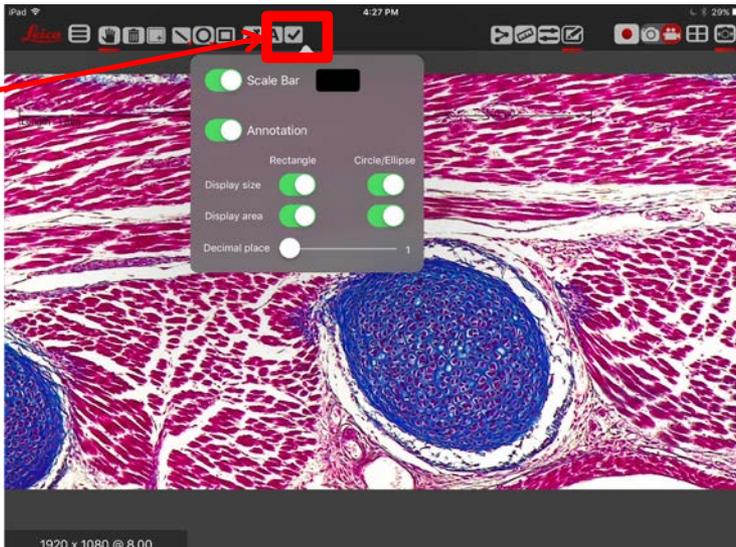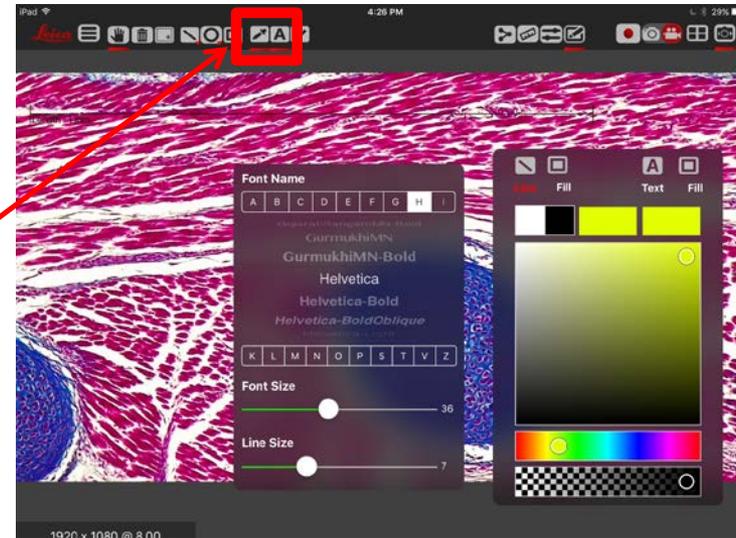

Color scheme

Font

On/Off

Still and Video Capture

- Select Still or Video Capturing





Acquire

Still or Video mode
(slider selection)



Share

- Share the image from Live Image View or send to device photo library or another collection.

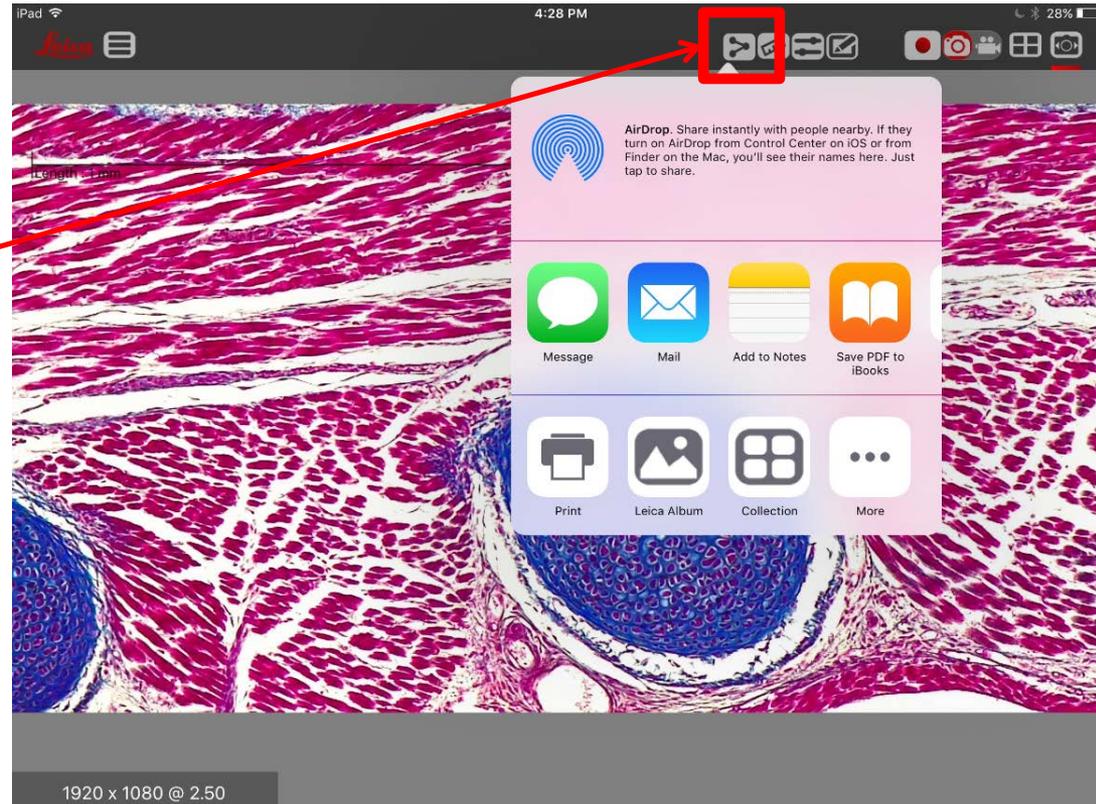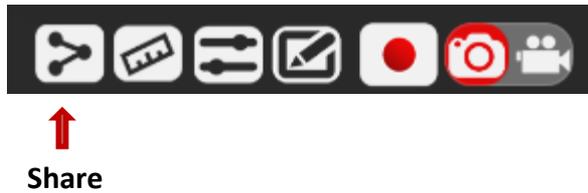


Gallery View

- Return to Gallery View



↑
Gallery View

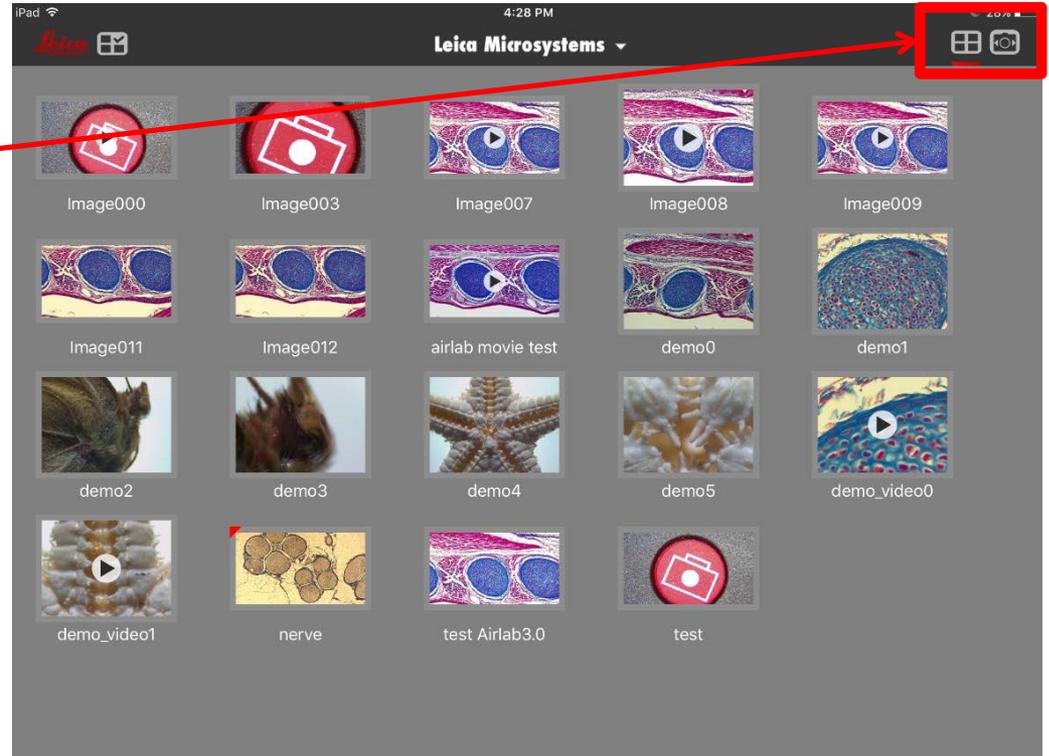

Select a Gallery Image

- Select a thumbnail image for an action




Select an image



Trash


Share

Select all or none

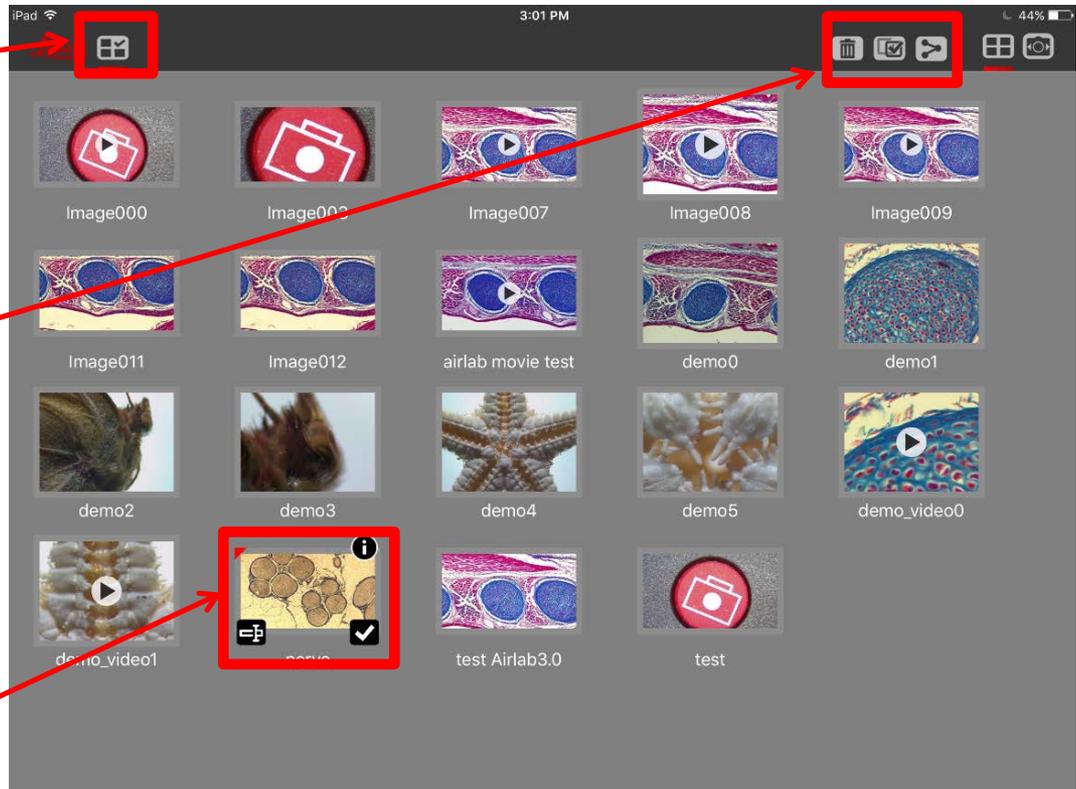

View Image Data

Indicates image contains annotations

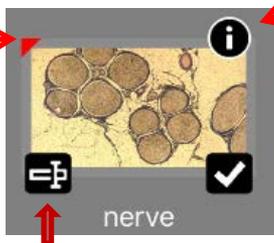

Change image name

Select a Gallery Image


- Tap a Gallery Image to view in full screen for Sharing, Image adjustments, or annotations



↑
Share

↑
Image adjustments

↑
Annotate

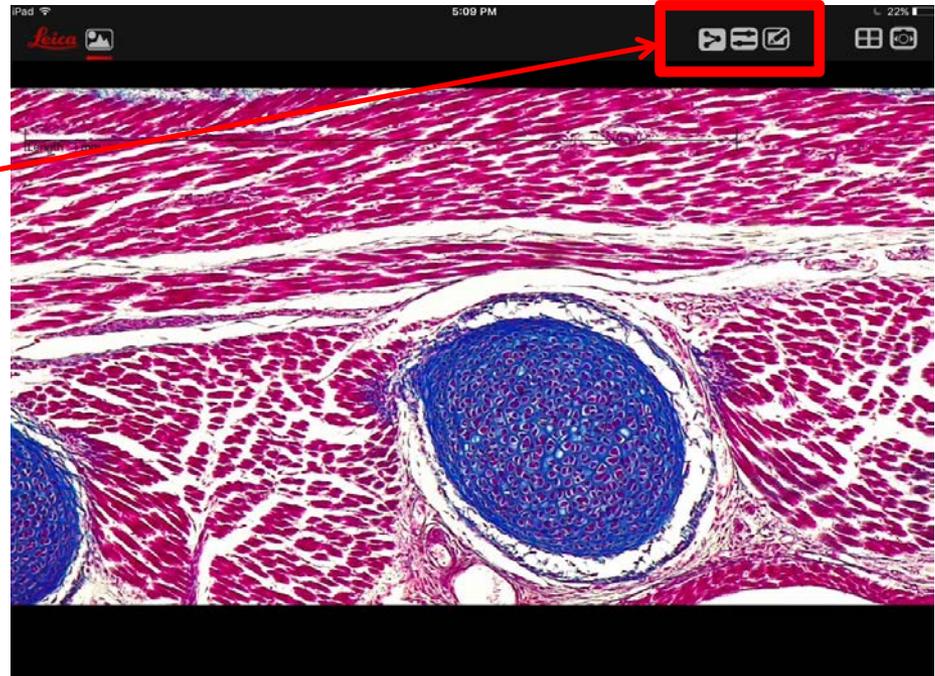

From Eye to Insight



Leica AirLab Icon Guide